# ADA USB/Ser

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### Allgemeines

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## Operating manual

### General information

The ADA USB/Ser is an adapter for the connection of WTW meters with a serial interface to the USB port of a Personal Computer. Thus you can operate meters from the PC (e. g. with the aid of the Windows program MultiLab pilot), transfer data from the meter to the PC, and update meter software.

The following meters are suitable:

- all handheld meters of the series 330/340, 330i/340i and 350i
- all portable meters of the series 197 and 197i
- all inoLab Level 2, Level 3 laboratory meters as well as inoLab series 730/735/750 and terminal 740
- handheld photometers pFotoFlex and pFotoFlex Turb (direct connection or via the LabStation)
- handheld turbidimeter Turb 430 IR (direct connection or via the LabStation)
- all inoLab Level 3 / series 740 active multifunction boxes.  
The multifunction boxes are supplied with power via the USB port. Therefore, no separate connection (e. g. via a Y cable or similar) is necessary.

### System requirements

- PC with Pentium processor or higher with at least one free USB port and a CD-ROM drive
- Windows 2000 / XP / Vista / 7.

### Installation

- Insert the enclosed installation CD, "Driver ADA USB/Ser", in the CD drive of your PC
- Install the USB driver on the PC.
- Follow the Windows installation instructions as necessary.
- Now you can connect the WTW meter to the 9-pin Sub-D socket of the ADA USB/Ser via the AK 340/B or AK 540/B cable.  
The LabStation can be connected using the zero modem cable (LabStation scope of delivery)
- The meter is listed as a virtual COM interface among the connections in the Windows instrument manager.

### Note:

Possible problems with older versions of the MultiLab pilot software can be eliminated by an update to the current version. This is available on the Internet under [www.wtw.com](http://www.wtw.com).

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